



CENTRIFUGE SAFETY



The centrifuge is a commonly used tool in laboratory research. It uses centrifugal force to separate substances in liquid or solid media according to particle size and density differences.

Hazards presented by all centrifuges, including microcentrifuges, if used and/or maintained improperly include:


- **Physical hazards:** Mechanical stress, metal fatigue, and corrosion of the rotor over time.
- **Exposure hazards:** Aerosolization of biological, chemical, or radioactive materials.

STANDARD OPERATING PROCEDURE GUIDANCE

The following information may be integrated into a lab-specific standard operating procedure (SOP) for centrifuge use.

1. PLANNING FOR USE

- Complete lab-specific training for the centrifuge.
- Wear appropriate PPE: Including safety eyewear, gloves, lab coat, and appropriate street clothing (i.e., long pants and closed-toe shoes). Ensure gloves are compatible with hazard(s).
- Inspect centrifuge:
 - Ensure tubes are rated for intended use (speed, temperature, and chemical resistance).
 - Rotor compatible with centrifuge and seated on drive correctly.
 - Rotor and safety cups/buckets free of cracks and deformities.
 - Rotor O-ring not cracked, missing, or worn.
 - Safety cups/buckets attached correctly and able to move freely.

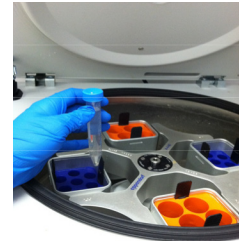

⚠ Contact a qualified technician if inspection identifies centrifuge components requiring repair or replacement.

- Prepare centrifuge tubes for loading:
 - Inspect centrifuge tubes before use.
 - Follow manufacturer's filling limits for tubes. Do not overfill or underfill tubes.
 - For biohazardous materials, disinfect outside of tubes prior to removal from biosafety cabinet and loading into rotor.

⚠ When centrifuging hazardous materials, use tightly capped tubes and/or sealable safety cups or rotors that can be loaded and unloaded in a fumehood or biosafety cabinet, depending on hazard.

- Use in-line filter for high speed centrifuges and ultracentrifuges to prevent contamination of vacuum pump and pump oil. Provide secondary containment for vacuum pump.

2. CENTRIFUGE OPERATION

- Balance Centrifuge: See the following link for tips on centrifugation without use of a balance tube.
<http://www.stanford.edu/dept/EHS/prod/researchlab/lab/centrifuge.html>
- Start run
 - Do not leave centrifuge unattended until full operating speed is reached and appears to be running safely without incident.
 - Stop centrifuge immediately if you notice any unusual noises or shaking. Confirm rotor is balanced.

⚠ To prevent rotor failure, do not exceed maximum speed and maximum mass limits for the rotor. You must reduce rotor speed if sample density calculations indicate maximum mass limits will be exceeded; contact manufacturer for guidance.

3. SAMPLE REMOVAL

- Stop run: Ensure centrifuge comes to complete stop before opening cover.


⚠ When centrifuging hazardous materials, wait at least 10 minutes after run to allow aerosols to settle before opening centrifuge.

- Check for leaks/spills: In samples, rotor, safety cups/buckets, and centrifuge well.
- Open sealable tubes/safety cups/rotors: Wear appropriate PPE and open inside fumehood or biosafety cabinet, depending on hazard.

CENTRIFUGE MAINTENANCE

1. PREVENTIVE MAINTENANCE

- Establish preventive maintenance schedule: Including regular cleaning of centrifuge interior to prevent damage and avoid costly repairs. Reference centrifuge operator's manual or contact manufacturer for guidance.

 Equipment repair and adjustments shall only be conducted by qualified service technicians.

- Maintain log book: For all high speed centrifuges and ultracentrifuges include run dates, durations, speeds, total rotor revolutions, and notes on rotor condition.

 Retire rotors after manufacturer's recommended life span except where annual stress test demonstrates absence of structural flaws. Note: Rotor life span may be reduced or warranty voided if autoclaved; contact manufacturer for guidance.



*Mechanical failure of an ultracentrifuge
By Permission of Cornell University*

2. CENTRIFUGE DISPOSAL

- For biohazardous materials, clean/disinfect centrifuge and remove/cross out biohazard sticker. Attach note on centrifuge describing what has been done.
- For radioactive materials, request radiation safety survey and signage before disposal of centrifuge. Contact EHS at (657) 278-7233.

EMERGENCY RESPONSE


1. MECHANICAL FAILURE

- Turn off centrifuge immediately and unplug power cord.

 Do not use centrifuge again until inspected by qualified service technician.

2. HAZARDOUS MATERIAL SPILL/EXPOSURE

- Turn off centrifuge immediately.

 Keep centrifuge cover closed for at least 30 minutes to reduce aerosolization of hazardous material.

- Follow spill, exposure, and incident reporting instructions.
 - http://ehs.fullerton.edu/_documents/emergencyinfo/chemicalspill.pdf
 - For radioactive materials: Call Environmental Health and Safety at (657) 278-7233 or (657) 278-2515.

REMEMBER:

- Conduct rotor or safety cup/bucket cleanup in nearest biosafety cabinet or fumehood, depending on hazard.
- Use tongs or forceps to avoid contact with sharp-edged debris.
- Avoid alkaline cleaners for aluminum centrifuge components.
- Avoid abrasive wire brushes for cleaning.

IMPORTANT CONTACTS

- Principal Investigator: _____(NAME), _____(PHONE).
- Equipment Service Technician (Chemistry/Biochemistry): Phil Berriman (657) 278-3662.
- Equipment Service Technician (Biology): Aaron Daveler (657) 278-2463.
- Environmental Health and Safety: (657) 278-7233.